How to ask ZMAN License Code

1. Download the latest version of ZMAN software and runtime engine at below links.

ZMAN software download

(path: www.zivelab.com - Support - Software download menu)

LabView runtime engine download

Cautions:

- 1) The license code is tied to a PC. Please perform the following steps on the PC which the ZMAN will be installed on.
- 2) The 32 bit LabView runtime engine 2013 is needed even if your OS is 64bit. The LabView runtime engine for 64 bit OS is not available for this software.

2. Run the ZMAN software.

	Open	News from WonATech
	Empty Project	 News from Zive Lab About ZMAN Request Supports Getting Started User Manual(2.3)
	🛅 Browse	
IVE ZMAN™ 2.3	Tools	Unknown
Impedance Spectroscopy nalysis & Presentation	za) Model Editor	Unknown Activation required

3. If you click 'Activate' button in red box, below window will pop out. Click 'Request Code' button.

Enter your license	e code,
If you want to ac	tivate the software, click the "Request Code
button. With the software without	license code, you can use all features of the restriction.
Places and that t	Alter - Ale
LICENCE PROFESSION FRANK	THE PERSON AND AND AND ADDRESS OF A DESCRIPTION AND ADDRESS ADDRES
you need a valid	license for each computer you wish use it.
you need a valid	license for each computer you wish use it.
viease note that t you need a valid License Code :	inis software is licensed per computer, so license for each computer you wish use it.

4. Click 'Select WonATech' button.

If you are using one of Zahner's instruments, click 'Select Zahner' button.

Please select your sales re	epresentative.
f you have Zahner produ	ucts, you may get a license code to
Zahner" button.	mer. In such a case, click the "Select
)therwise, please select V	WonATech.

5. Fill out the below windows and 'OK' button.

	* Required
*First Name	*Last Name
*E-Mail Address	
*Phone Number	Fax Number
*Company/Institute	Department
*Street Address 1	
Street Address 2	
*Country	*Zip/Postal Code
Note	
Your right to privacy is a need for reasonable cont with us. Therefore, this st guideline, and is not limit not sell your information goal in gathering informa our relationship with you	priority to us. We recognize your rol of information that you share atement is a corporate-wide ted to our online business. We do to outside organizations. Our ation is to consistently improve , our customer.

6. Click 'Send Email', then we will receive your e-mail asking a ZMAN license code.

Or, if your computer is not connected to the internet, please copy this information and sent it to sales@wonatech.com.

User Info	×
Please e-mail the followings to your sales re (sales@wonatech.com). You will receive an license code from the representative. You can also find it in the clipboard. Please any text editor and e-mail it later. Product: ZMAN2x	epresentative e-mail with a valid paste(CTRL-V) it in
Host Name: DIANE PG- Host ID: BOTA 6653 0000 077A	
First Name: 1 Last Name: 1 E-Mail Address: 1@1 Phone Number: 1 Fax Number: Company/Institute: 1 Department Street Address1: 1 Street Address2: Zip/Postal Code: 1 Country: 1 Serial Number:	
Before	Send Email

7. After you receive a license code, go to step 3 to enter the license code you received.

Enter your licen	se code.
If you want to a button. With th software withou Please note that	activate the software, click the "Request Code e license code, you can use all features of the it restriction. t this software is licensed per computer, so
you need a vali	a license for each computer you wish use it.
License Code :	3144-5618-EECF-398D

8. Now your ZMAN is activated.

	Open	News from Won ATach
4	Empty Project	 Discussion Forums Knowledge Base Request Supports Getting Started User Manual(2:2 F3)
	🛅 Browse	
ZMAN TM 2.2	Tools	test1 test2
F3 patch Impedance Spectrescopy Analysis & Presentation	25a Model Editor	test company 3144-5618-EECF-398D